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Instruction manual for a hydrometer for measuring alcohol in volume percentage (% vol. Gay-Lussac) 20° C

The alcoholmeter for % vol. is for the determination of the alcoholic content in water. The instrument indicates therefore only a pure alcohol- water mixture. It cannot be used for measuring liqueurs of fluids in which sugars or other materials besides water and alcohol. These would lead for the falsification of the advertisement.

Since the alcoholmeter is fragile, it should be treated carefully, dry and clean. The jar from glass should in the diameter at least 1,5 – 2,0 cm broader is than the diameter of the hydrometer. It must be likewise clean and dry and it must stand vertically. The testing liquid (alcohol water mixture) should be as clean as possible all other parts should be removed previously with sieves of filters.

The Alcoholmeter is touched at the thin end above the **incasuring**-scale carefully and slowly put into the liquid. It may immerse thereby at most 5 mm more deeply than the natural swimming position. Please give your attention to the distance between Alcoholmeter and jar. An adhesion of touch is to avoid. After immobilization of the Alcoholmeter (30sec) the results can be read.

Please bring the eye densely underneath the meniscus so that the stem becomes an elliptical plain around it. After lifting of the eye slowly becomes these to a line. Where this line meets the scale you found the reading position like enclosed drawing.